

COMPOUNDING MADE SIMPLE

UNGUATOR E/S

Streamline compounding in your practice with the help of the Unguator E/S.

This electronic mortar and pestle offers quick and consistent preparations while maintaining quality and minimizing contamination. This unit allows the operator to simply yet effectively mix topical compounds with the push of a button.

Simply weigh your ingredients into the Unguator Jar and mount it to the machine to be mixed. Next select your mix duration and your speed settings then press start. The automated mix will now begin.

While the mix is running you can take this time to do other valuable tasks within your pharmacy.

Once the mix is complete, remove the jar from the machine; put your prescription label on the jar, and give the prescription to the patient. It's that simple!



Electronic Mortar and Pestle!

www.hescorp.net | (405) 329-6810